# PolyBranch HCR™ Rapid RNA FISH Kit(Plant)

**Catalog Number:HKR11Z**

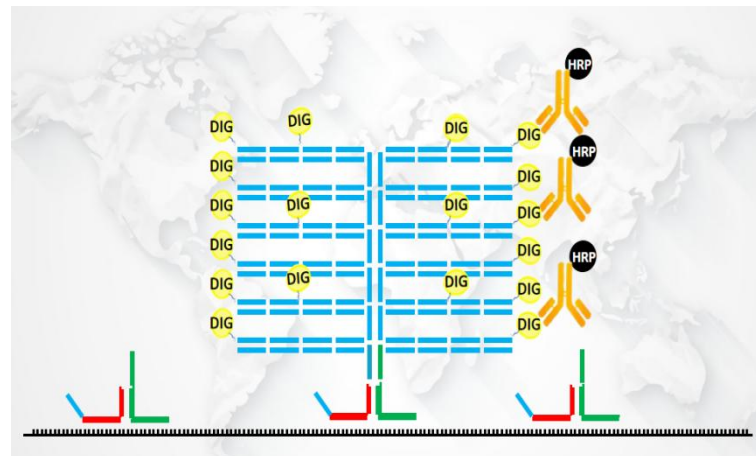
## 【Product Information】

| Product Name                                             | Catalog No | 20T   | Storage                         | Shelf Life |
|----------------------------------------------------------|------------|-------|---------------------------------|------------|
| Wash Buffer (powder, dissolve in 10L ddH <sub>2</sub> O) | HKR11-1    | 10L   | Short-term:4℃<br>Long-term:-20℃ | 12 months  |
| Proteinase K (100x)                                      | HKR11-2    | 20uL  |                                 |            |
| Pectinase/Cellulase Mixture                              | HKR11-3    | 2mL   |                                 |            |
| AP-Mouse Anti-Digoxin (100x)                             | HKR11-4    | 20uL  |                                 |            |
| NBT Chromogen Solution A (20x)                           | HKR11-5    | 100uL |                                 |            |
| NBT Chromogen Solution B (20x)                           | HKR11-6    | 100uL |                                 |            |
| Blocking Buffer                                          | HKR11-7    | 2mL   |                                 |            |
| Probe:                                                   | HKR11-8    | 2mL   |                                 |            |

## 【Product Description】

RNA rapid hybridization kit, suitable for in situ hybridization of RNA larger than 300bp, with simple and convenient operation and short experimental time. The reagent kit contains multiple pairs of trigger probes. When two probes bind to adjacent positions of the target sequence, one probe triggers the other probe to expose the amplified sequence. The amplified sequence triggers a branched HCR to form a large molecular nucleic acid aggregate, and each nucleic acid strand is labeled with digoxin. The antibody recognizes digoxin for color

development.

## 【Probe Information】

## 【Tissue Fixation】

| Tissue Type | Animal                            | Plant                                              | Frozen Samples                                                                        | Cell Climb Slides  | Cells                                                                        |
|-------------|-----------------------------------|----------------------------------------------------|---------------------------------------------------------------------------------------|--------------------|------------------------------------------------------------------------------|
| Treatment   | Fix at RT for 12h, paraffin embed | Vacuum fix for 1h, RT fix for 12h, paraffin embed. | Dehydrate in 15% sucrose at 4°C for 8h, then in 30% sucrose at 4°C for 8h, OCT embed. | Fix at 4°C for 2h. | scrape off cells, fix in 4% PFA at 4°C for 2h, wash with PBS, agarose embed. |
| Type        | mRNA                              | lncRNA                                             | circRNA                                                                               | miRNA              | rRNA                                                                         |
| Treatment   | Fix at RT for 12h                 | Fix at RT for 12h (<300bp: 24h)                    | Fix at RT for 12h                                                                     | Fix at RT for 12h  | Fix at RT for 12h (<300bp: 24h)                                              |

## 【Storage and Shipping】

Ship on wet ice; store at -20 °C for long-term or at 4 °C for short-term use. Shelf life: 6 months.

## 【Protocol】

### 1. Deparaffinization and Rehydration

Immerse slides sequentially in: Xylene I (15 min) → Xylene II (15 min) → Xylene III (15 min) → 100% ethanol (10 min) → 90% ethanol (10 min) → 80% ethanol (10 min) → 70% ethanol (10 min) → Rinse with distilled water.

### 2. Enzyme Repair

After slides are completely dry, draw a hydrophobic circle around the tissue using a histology pen (recommended: HKR14P In Situ Hybridization Pen). Place slides horizontally in a hybridization oven or humidified chamber. Add 100 µL of Pectinase/Cellulase Mixture onto the tissue and incubate at 37 °C for 30 min. Rinse with distilled water to stop the reaction. Remove excess liquid, then add 100 µL of Proteinase K repair solution (1X) and incubate at 37 °C for 30 min. Rinse with distilled water to stop the reaction.

### 3. Blocking

Remove excess liquid from slides. Add 100 µL of Blocking Buffer per slide and incubate at 37 °C for 30 min. Wash once for 5 min. **Wash steps: Place slide racks in a wash tank, add wash buffer (ensure samples are submerged), and shake at 60 rpm for 5 min.**

### 4. Probe Hybridization

Remove excess liquid from slides. Add 100 µL of probe per slide and incubate at 37 °C for 2 – 3 h (maintain humidity to prevent drying). Wash 5 times, 5 min each. **Wash steps: As described above.**

### 5. AP-Mouse Anti-Digoxin

Remove excess liquid from slides. Add 100 µL of AP-Mouse Anti-Digoxin (1X) per

slide and incubate at 37 °C in a humidified chamber for 40 min. Wash 5 times, 5 min each. **Wash steps: As described above.**

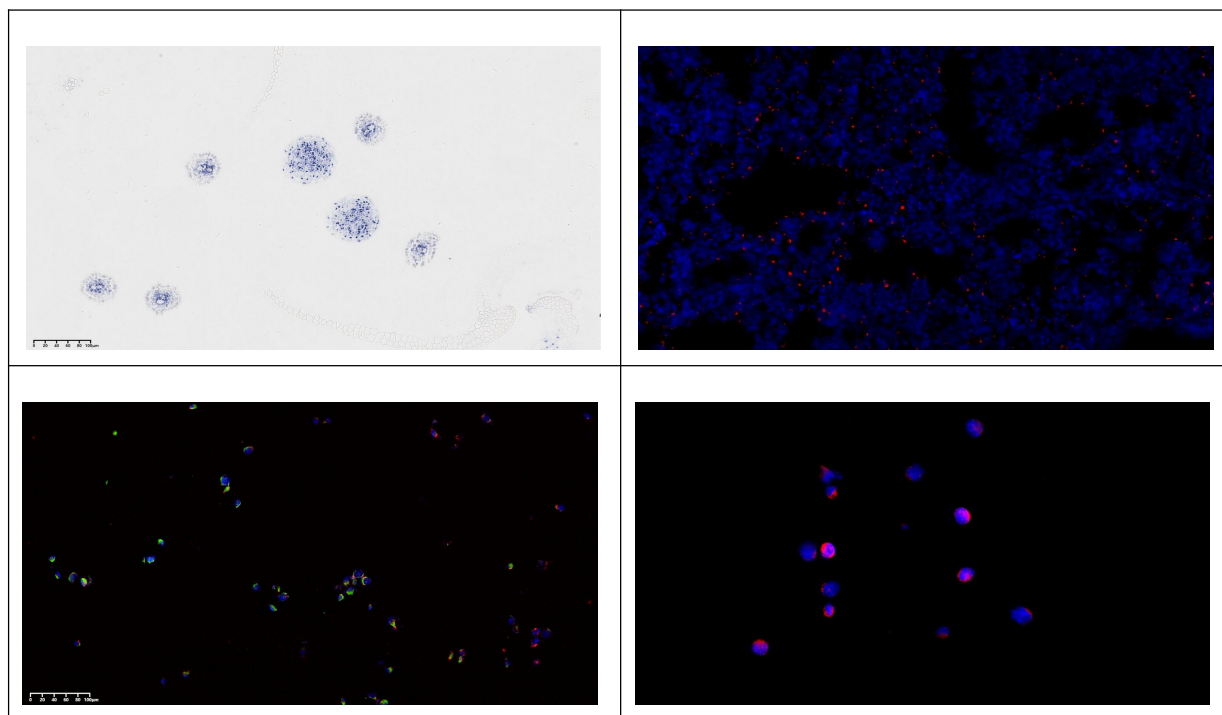
## 6. Chromogenic Reaction

Prepare the chromogenic solution by adding 50  $\mu$ L of NBT Chromogen Solution A and 50  $\mu$ L of NBT Chromogen Solution B to 1 mL of deionized water. Add 100  $\mu$ L of the mixture per slide and incubate at RT for 30 min. Rinse with distilled water to stop the reaction. Air-dry slides and mount with mounting medium.

### 【Precautions】

1. For research use only.
2. Wear lab coats and disposable gloves for safety.

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